

Membrane Degasification Chemical-free alternative for the pharmaceutical industry

Technology for the Removal of Dissolved Gases

In treating water for the production of pharmaceuticals, the concentration of gases dissolved in the water plays an important role. In many cases the CO₂ content in the raw water has to be further reduced to obtain purified water with a conductivity of < 1.3 μS/cm (at a temperature of 25°C).



Contactors in an 8m³/h ROCEDIS plant

Historically, NaOH dosing has been used to control CO₂. With chemical dosing CO₂ is converted into a carbonate, which can be removed by the RO. The latest state of the art technology for this task is membrane degasification where chemicals are not needed.

Design / Function

ONDEO Industrial Solutions - Hager + Elsässer, is an OEM who now specifies Liqui-Cel Membrane Contactors for CO₂ removal for pharmaceutical water systems. These system designs place the

membrane degasification system downstream of the RO system and before an EDI system. See the flow diagram on page 2.

A Polypropylene Hollow Fiber membrane is used to put a gas and liquid phase in direct contact with each other. Because the fiber is hydrophobic, the liquid will not penetrate the pore. A strip gas or vacuum used on the inside of the hollow fiber lowers the partial pressure of the gas phase


which causes the gases to diffuse from the liquid phase through the membrane wall into the gas phase. The removed CO₂ is continuously swept out of the contactor by the air sweep.


Hollow fiber membrane contactors maximize the surface contact area for efficient, chemical-free, CO₂ removal.

Advantages

- Chemical free
"no added substances"
- Continuous operation
- Compact design
- Easy to expand due to modular design
- Reduction in operating costs

Since 1996 Hager + Elsässer has been integrating Membrane Degasification in their pharmaceutical water treatment systems. World-wide more than 90,000 m³/h (400,000 gpm) of water have been treated with Membrane Contactors.

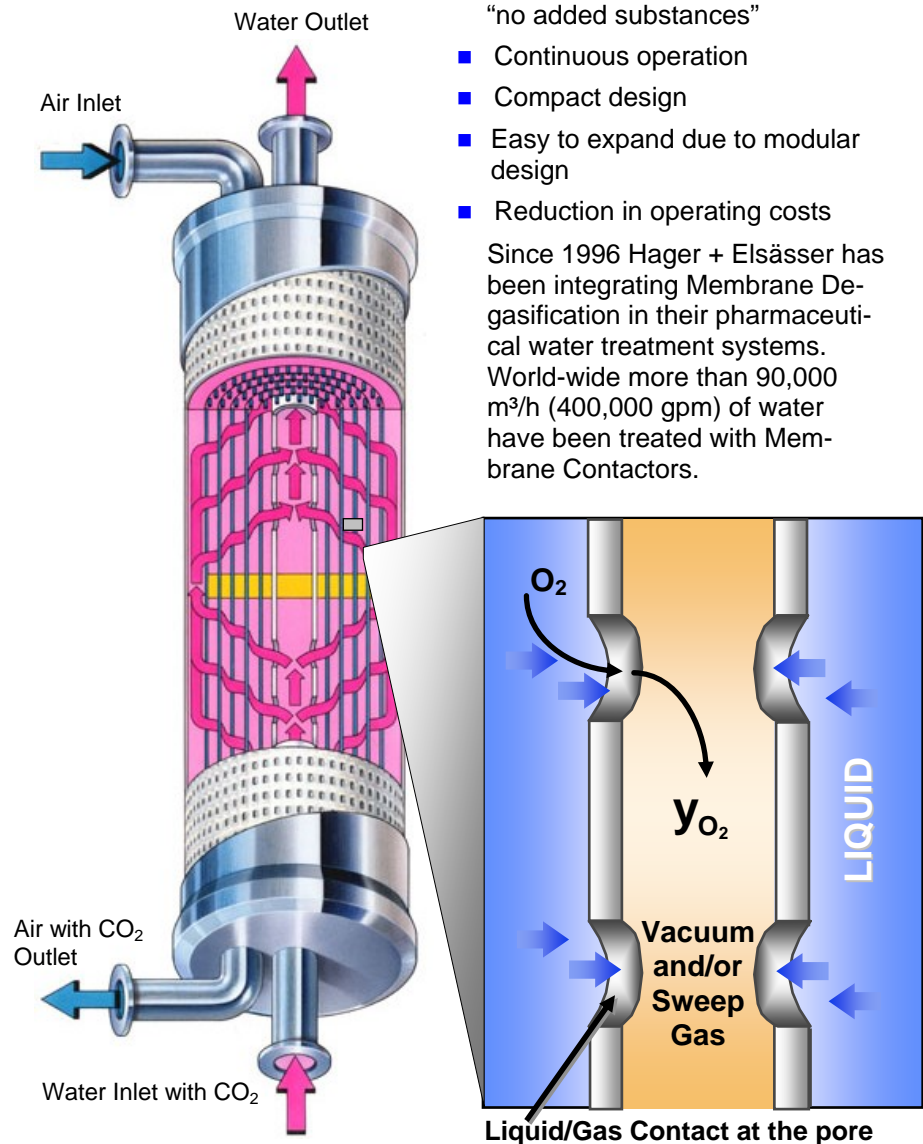

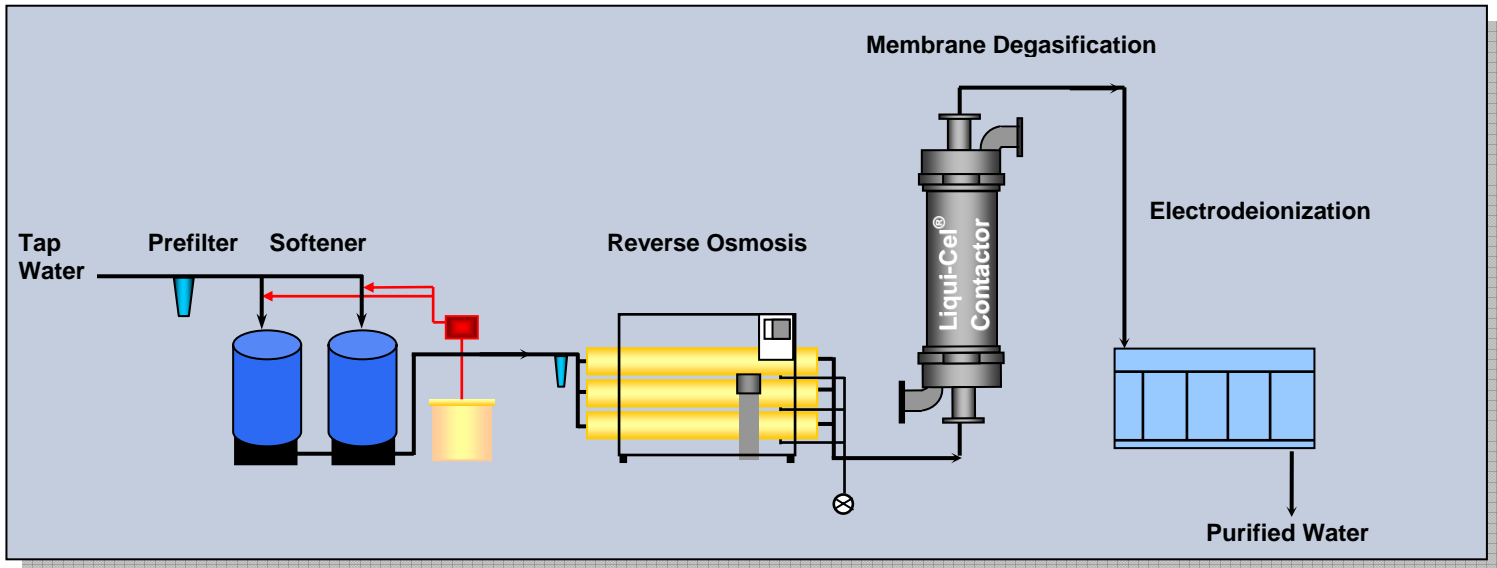

Flow diagram of the plant design to produce Purified Pharmaceutical Grade Water per US and European Pharmacopoeia Requirements

Produced in conjunction with ONDEO Industrial Solutions GmbH - Hager+Elsässer Centre of Excellence. For more information, visit www.hager-elsaesser.com

This product is to be used only by persons familiar with its use. It must be maintained within the stated limitations. All sales are subject to Seller's terms and conditions. Purchaser assumes all responsibility for the suitability and fitness for use as well as for the protection of the environment and for health and safety involving this product. Seller reserves the right to modify this document without prior notice. Check with your representative to verify the latest update. To the best of our knowledge the information contained herein is accurate. However, neither Seller nor any of its affiliates assumes any liability whatsoever for the accuracy or completeness of the information contained herein. Final determination of the suitability of any material and whether there is any infringement of patents, trademarks, or copyrights is the sole responsibility of the user. Users of any substance should satisfy themselves by independent investigation that the material can be used safely. We may have described certain hazards, but we cannot guarantee that these are the only hazards that exist.

Liqui-Cel, Celgard, SuperPhobic and MiniModule are registered trademarks and NB is a trademark of Membrana-Charlotte, A division of Celgard, LLC and nothing herein shall be construed as a recommendation or license to use any information that conflicts with any patent, trademark or copyright of Seller or others.

©2007 Membrana – Charlotte A Division of Celgard, LLC (TB38 Rev1 10/05)

Membrana – Charlotte
A Division of Celgard, LLC
13800 South Lakes Drive
Charlotte, North Carolina 28273
USA
Phone: (704) 587 8888
Fax: (704) 587 8585

Membrana GmbH
Oehder Strasse 28
42289 Wuppertal
Germany
Phone: +49 202 6099 - 658
Phone: +49 6126 2260 - 41
Fax: +49 202 6099 -750

Japan Office
Shinjuku Mitsui Building, 27F
1-1, Nishishinjuku 2-chome
Shinjuku-ku, Tokyo 163-0427
Japan
Phone: 81 3 5324 3361
Fax: 81 3 5324 3369

MEMBRANA
Underlining Performance
www.membrana.com
www.liqui-cel.com

A **POLYPOR** Company